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H_IL-5 Reporter 293 Cell Line

H_IL-5 Reporter 293 细胞系

For research use only!

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Reagent	Specification	Manufacturer/Catalogue No.
Puromycin	25 mg	Genomeditech/GM-040401-1
Blasticidin	10 mg	Genomeditech/GM-040404-1
Hygromycin	1 g	Genomeditech/GM-040403-1
G418	10 mg	Genomeditech/GM-040402-1
Pen/Strep	100 mL	Thermo/15140-122
Fetal Bovine Serum	500 mL	ExCell/FSP500
DMEM	500 mL	gibco/C11995500BT
96 well round cell culture plate	Plate	PakGent/CL-PT096
96-well White Opaque Plate	Plate	Thermo/236108
96-well U-bottom Plate	Plate	Saining/1014010
Human IL-5 Protein; His Tag	/	Genomeditech/GM-88539RP
Recombinant Human IL-5 Protein	5 µg	Sinobiological/15673-HNCE
Mepolizumab (anti-IL-5)	100 µg	Aladdin/Ab170275-100µg
Anti-IL-5R hIgG1 Antibody(Benralizumab)	/	Genomeditech/GM-46851AB
Anti-IL-3R Beta/IL3RB Antibody (APC), Mouse Monoclonal	25 Test	Sino Biological/10516-MM01-A
GMOne-Step 2.0 Luciferase Reporter Gene Assay Kit	/	Genomeditech/GM-040513C

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Equipment	Manufacturer/Catalogue No.
细胞计数仪	ThermoFisher Scientific/Countess 3
酶标仪	Moleculardevices/SpectraMax L

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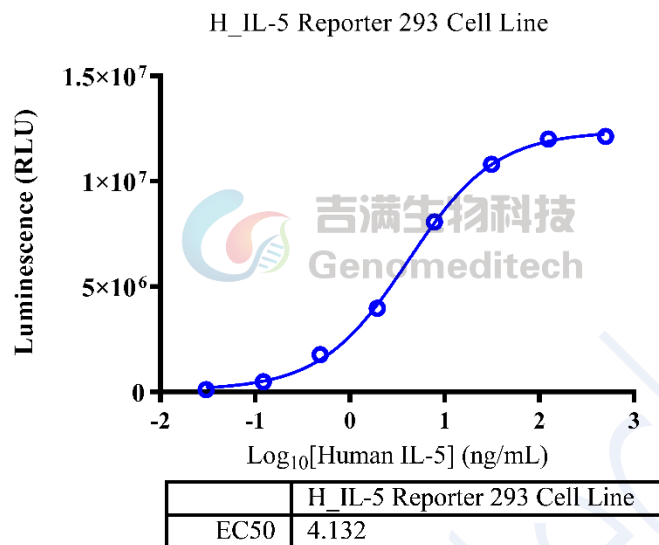
Fig 2. Response to Human IL-5 Protein; His Tag. The H_IL-5 Reporter 293 Cell Line (Cat. GM-C06723) at a concentration of 1.5E4 cells/well (96-well format) was stimulated with serial dilutions of Human IL-5 Protein; His Tag (Cat. GM-88539RP) in assay buffer (DMEM + 1% FBS + 1% P.S) for 16 hours. The firefly luciferase activity was measured using the Luciferase Reporter Assay Kit (Genomeditech). The maximum induction fold was approximately [105]. Data are shown by drug mass concentration.

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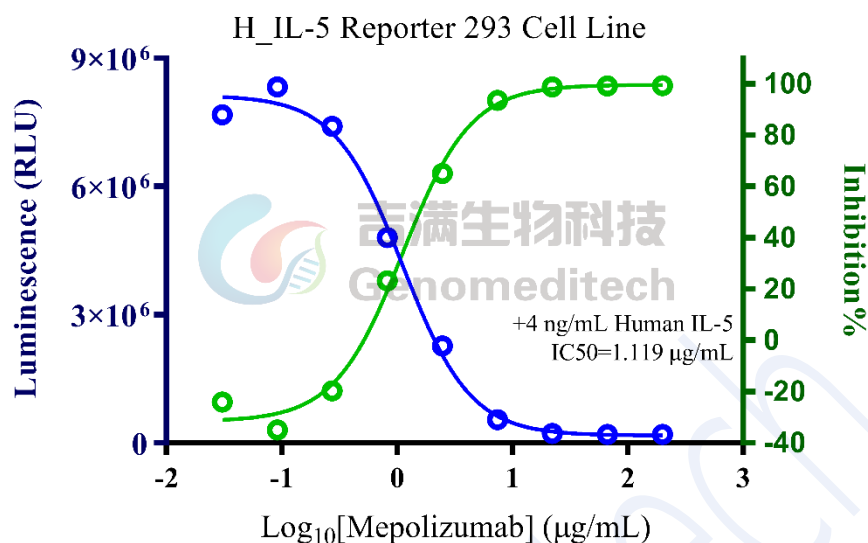
Fig 3. Response to Mepolizumab (anti-IL-5). Serial dilutions of Mepolizumab (anti-IL-5) (Aladdin/Ab170275-100μg) was incubated with 0.4 ng/well of Human IL-5 Protein; His Tag (Cat. GM-88539RP) for 1 hour in assay buffer (DMEM + 1% FBS + 1% P.S). After pre-incubation, add the mixture to the H_IL-5 Reporter 293 Cell Line (Cat. GM-C06723) at a density of 1.5E4 cells/well in a 96-well format, and incubate for 16 hours. Firefly luciferase activity was measured using the Luciferase Reporter Assay Kit (Genomeditech) (left Y-axis, relative luminescence units, RLU), with inhibition percentages shown on the right Y-axis. Data are shown by drug mass concentration.

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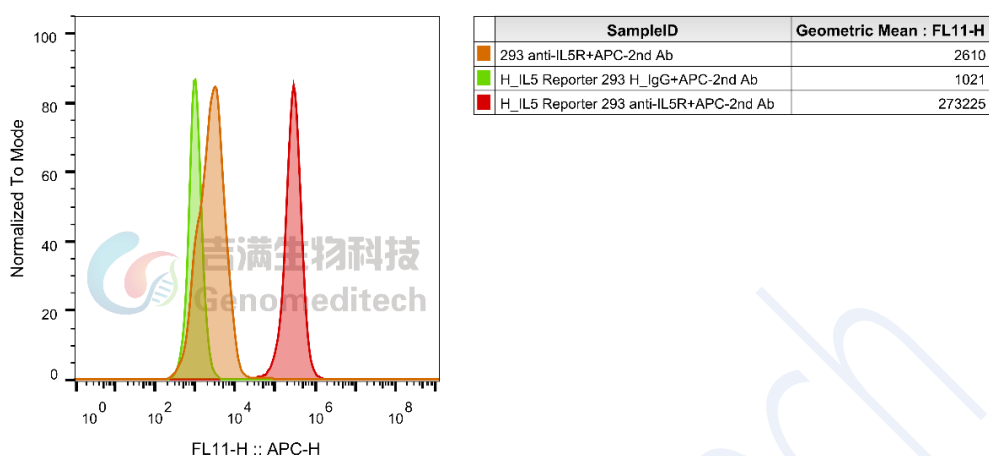
Fig 4. H_IL-5 Reporter 293 Cell Line (Cat. GM-C06723) was determined by flow cytometry using Anti-IL-5R hIgG1 Antibody(Benralizumab) (Cat. GM-46851AB).

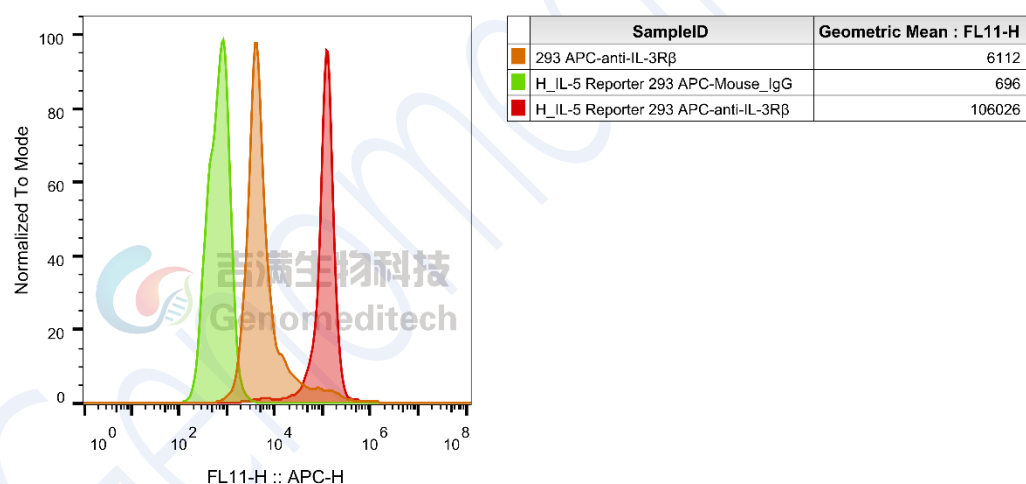
Fig 5. H_IL-5 Reporter 293 Cell Line (Cat. GM-C06723) was determined by flow cytometry using Anti-IL-3R Beta/IL3RB Antibody (APC) (SinoBiological/10516-MM01-A).

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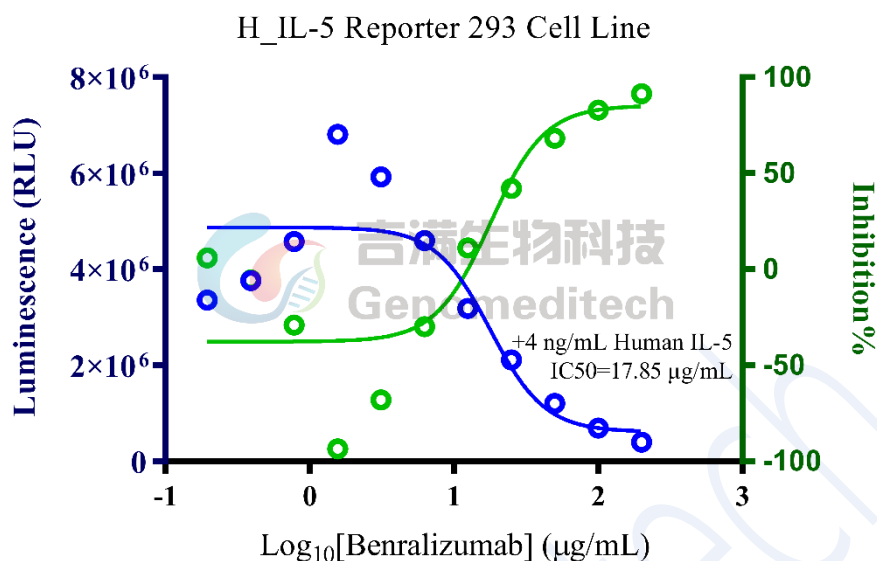
Fig 6. Response to Benralizumab. Serial dilutions of Anti-IL-5R hIgG1 Antibody(Benralizumab) (Cat. GM-46851AB) and 0.4 ng/well of Human IL-5 Protein; His Tag (Cat. GM-88539RP) were added to the H_IL-5 Reporter 293 Cell Line (Cat. GM-C06723) at a density of 1.5E4 cells/well in a 96-well format, and incubated for 16 hours. Firefly luciferase activity was measured using the Luciferase Reporter Assay Kit (Genomeditech) (left Y-axis, relative luminescence units, RLU), with inhibition percentages shown on the right Y-axis. Data are shown by drug mass concentration. Notably, low concentrations of the anti-IL-5RA antibody may exhibit an activation effect.

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Zabeau L, Van der Heyden J, Broekaert D, Verhee A, Vandekerckhove J, Wu SJ, Chaiken I, Heinrich P, Behrmann I, Tavernier J. Neutralizing monoclonal antibodies can potentiate IL-5 signaling. *Eur J Immunol.* 2001 Apr;31(4):1087-97. doi: 10.1002/1521-4141(200104)31:4<1087::aid-immu1087>3.0.co;2-q. PMID: 11298333IF: 4.1 Q2 .

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IL-4/IL-13	
IL-4 Reporter Cell Line	IL-4/IL-13 Reporter 293 Cell Line
IL-4/IL-13 Reporter 293 DDX35TM Cell Line	Cynomolgus_IL4R CHO-K1 Cell Line
H_IL4R CHO-K1 Cell Line	
Anti-IL-4R hIgG1 Antibody(12B5)	Anti-IL4R hIgG4 Antibody(Dupilumab)
Anti-IL4R hIgG4 Reference Antibody (Dupbio)	
Cynomolgus IL-4R alpha Protein; His Tag	Human IL-4R alpha Protein; His Tag
Human IL-4R alpha Protein; mFc Tag	
TSLP:TSLPR	
H_TSLP Reporter Cell Line	H_TSLPR CHO-K1 Cell Line
Anti-H_TSLPR hIgG1 Antibody	Anti-TSLP hIgG2 Reference Antibody(Tezbio)
Anti-TSLP hIgG2 Antibody(Tezepeumab)	
Cynomolgus TSLP Protein; His Tag	Human TSLP Protein; His Tag
IL-5	
H_IL-5RA CHO-K1 Cell Line	H_IL-5RA HEK-293 Cell Line
Anti-IL5 hIgG4 Antibody(Reslizumab)	Anti-IL-5R hIgG1 Antibody(Benralizumab)

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